

STANDARD OPERATING PROCEDURES
DIVISION OF COMPARATIVE MEDICINE
UNIVERSITY OF SOUTH FLORIDA

SOP#: 1166

Date Issued: New

Date Revised: 7/19

Page 1 of 2

TITLE: Sentry Digital Ultrasonic Cleaner

SCOPE: Animal Care Personnel

heavy bioburden is present it is recommended to use a detergent safe cleaning agent. If a heavy bioburden is present it is recommended to use a detergent safe cleaning agent.

2. Open the lid and place items in the tank. Add water to the tank to cover the surface of the items. The water level should stay between the MIN and MAX mark.
3. Cleaning agents may be added to the water to enhance cleaning.
4. Plug unit into electrical socket and observe green LED light on right side come on. PRESS the "ON" button to turn unit on. The digital display will show 15:0 (default cycle time is 15 min). Press the "ON" button again to start the cleaning cycle. To select a different cleaning cycle press "SET" and select from the preset cycles.
5. When the timer counts down to "000" the unit will turn off. The unit can be stopped/paused at any time by pressing the "OFF" button. Pressing the "ON" button will resume the cleaning and countdown.
6. Pressing the "TC" button turns on the heater function and illuminates the display.

g cycle

g the "OFF" button for more than 2 seconds turns the display off and puts the unit in standby mode.

